# SNAIGĖ

---

---

---

---

**MD14DC**

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Naudojimo instrukcija

**MEDICAL REFRIGERATOR**

Instruction for use

## GERBIAMIEJI!

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**DEAR CUSTOMER!**

CAREFULLY READ THE INSTRUCTION MANUAL. If instructions are not followed, there is a risk of injury, damage of the appliance and loss of the right to free warranty service.

The Manufacturer shall not be held liable for damage arising from the failure to observe the instructions contained in this manual.

Keep this manual for future reference, or to pass it over to the next user.

This instruction applies to more than one model and may contain inconsistencies.

**General information:**

The refrigerating device is a medicine refrigerator designed for professional storage of medicines and pharmaceutical preparations at a temperature of +5 °C.

The refrigerator is suitable for medicines, etc. for pharmaceutical preparations that need to be cooled and stored at a temperature of +2 °C to +8 °C.

The temperature-energy parameters of the refrigerating device meet the requirements of DIN 13277 (DIN 58245).

**Do not use refrigeration inside the appliance:**

- For cooling and storage of chemically unstable, flammable or alkaline substances.
- For cooling and storage of blood, plasma or other infused body fluids.

**GUIDELINES CONCERNING THE SAFETY OF USE****General safety requirements:**

- WARNING: Do not cover the ventilation openings of the appliance.
- WARNING: Do not use electrical appliances inside the storage compartments of the appliance, unless they are of the type recommended by the manufacturer.
- WARNING: Do not use any mechanical tools or other/chemical agents other than those recommended by the manufacturer to speed up the defrosting process.
- WARNING: When positioning the appliance, ensure the supply cord is not trapped or damaged.
- WARNING: Do not locate multiple portable socket-outlets or portable power supplies at the rear of the appliance.
- WARNING: Disposal of the appliance should be according to national rules.

WARNING: Do not damage the appliance refrigeration system. It contains the refrigerant gas R600a. If the refrigeration system is damaged:

- Do not use any open flame.
- Avoid sparks - do not turn on/off any electrical appliances or lighting fixtures.
- Immediately ventilate the room: air the room in which the appliance is placed for a few minutes (the size of the room for a product containing isobutane/ R600a must be at least 4 m<sup>3</sup>).

C-Pentane is used as blowing agent in the insulation foam and it is a flammable gas.

- This appliance is designed only for cooling and storing of pharmaceuticals.
- This appliance is not designed for storage of explosives, such as aerosols with flammable gases.
- It is prohibited to store petrol and other flammable liquids near the appliance.
- This appliance can be used by children aged from 8 years and above and persons with reduced physical, sensory or mental capabilities or lack of experience and knowledge if they have been given supervision or instruction concerning use of the appliance in a safe way and understand the hazards involved.

### **Requirements for the connection to the mains:**

- The appliance should ONLY BE CONNECTED TO AN EARTHED SOCKET INSTALLED IN ACCORDANCE WITH THE REGULATIONS. Make sure that the supply voltage corresponds with the voltage marked on the rating label. This product complies with all binding CE labelling directives. The earthed electrical socket by which the appliance is connected to the mains should be in an accessible place. It is a legal requirement that the appliance is properly earthed. The manufacturer will not be held liable for any damage or injury which may result from the failure to fulfil this requirement.
- It is mandatory to replace a damaged power cord with a proper one of the same type as the one used by the refrigeration appliance manufacturer to prevent damage to health and property. To avoid the exposure to danger, always have faulty cable be replaced only by the manufacturer, by our customer service or by a qualified person and with a cable of the same type.
- It is forbidden to readjust or modify any parts of the unit. It is crucial not to damage the capillary tube visible in the compressor recess. The tube may not be bent, straightened or wound. If the capillary tube is damaged by the user the guarantee will be void.
- Do not use adapters, multiple sockets and two-wire extension leads. If it is necessary to use the extension lead, it must be equipped with a protection ring and a single socket and must have a VDE/GS safety certificate. If an extension lead is used (with a protection ring and safety certificate), its socket must be located at a safe distance, away from the sinks, and must not be in a place where it could be flooded by water or waste water.
- The frequency and power of the electricity supply in your house must conform to the general data parameters of the appliance as it is shown in the product label.
- When positioning the appliance, be careful that the electrical cord isn't squeezed in order to avoid damaging it. Do not store heavy objects such as cooling devices or other devices next to the appliance in such a way that they could squeeze and damage the electrical cord. This can cause a short circuit and a fire.
- Make sure that the plug of the electrical cord is not squeezed by the back wall of the appliance or otherwise damaged. **A DAMAGED PLUG CAN BE THE CAUSE OF A FIRE!**

– **CAUTION!** If the installed interior lighting (LED) lamp is not functioning, contact the maintenance and repair representative for replacements. LAMPS CAN BE REPLACED ONLY BY THE MAINTENANCE AND REPAIR REPRESENTATIVE

 DO NOT TRY REPAIR THE APPLIANCE ON YOUR OWN. ALWAYS CONTACT AN AUTHORIZED TECHNICIAN. To avoid the exposure to danger, always have faulty cable be replaced only by the manufacturer, by our customer service or by qualified person with a cable of same type.

### **Requirements for safe use:**

**WARNING:** the device may only be operated with original accessories from the manufacturer or accessories from other suppliers approved by the manufacturer. Use of non-approved accessories is at the user's risk.

- It is forbidden to use a technically damaged appliance.
- If the appliance is unplugged (for cleaning, moving to another place, etc.), it may be repeatedly switched on after **15 min.**
- When performing activities, such as cleaning, maintenance or moving, the appliance must be completely disconnected from power supply (by pulling the plug out from the socket). **Do not pull on the cord, but hold the body of the plug instead.**
- Do not place heavy objects on the top of the refrigerating appliance.
- Do not place any switched on electrical devices on top of the appliance because this may cause ignition of plastic parts.
- Do not place any dishes with liquids on top of the appliance and do not keep flowers in vases or other liquid-filled vessels on the appliance.
- Do not climb on or sit on the appliance, do not lean on or hang on the appliance doors and do not allow children to do this.
- Do not use water spray and steam to clean the appliance.

 **The appliance should be transported in the vertical position;** do not tilt any more than 30°. THE MANUFACTURER WILL NOT BE RESPONSIBLE FOR ANY DAMAGE OF THE APPLIANCE THAT RESULTS FROM NON-COMPLIANCE WITH THE INSTRUCTIONS FOR TRANSPORTATION.

## **INSTALLATION AND OPERATING CONDITIONS**

### **Positioning:**


- Place the device in a dry, well-ventilated room with a minimum ambient temperature of +10 °C and a maximum of +35 °C.

**CAUTION!** Do not install this appliance in areas that are too humid or too cold. Place the appliance away from heat sources such as kitchen stove/oven, radiators, or direct sunlight. The appliance must not touch any pipes for heating, gas or water supply, or any other electrical devices.

- Do not cover the ventilation holes at the top of the appliances – it must be a good air circulation around the appliance. There should be a gap of at least 10 cm between the top of the appliance body and any furniture that may be above it. If this requirement is not followed, the appliance consumes more electrical energy and its compressor may overheat.

If the appliance is placed in a corner, a gap of at least 5 cm must be left between the appliance body and the wall.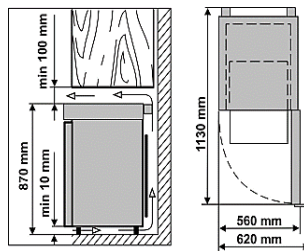

- The appliance must stand on a level surface and must not touch the wall. If necessary, regulate the height of the appliance by adjusting the levelling feet: by turning them clockwise – the front of the appliance rises, by turning them counter clockwise – it comes down. If the appliance is tilted slightly backward – the doors will close by themselves.

### Preparing for operation:

- It is recommended to prepare the appliance for operation with a helper.
- Remove package. Lift the appliance away from foamed polystyrene base. Therefore take out adhesive tapes from sides and simply lift the appliance up and take out the board. Strip adhesive tapes off and open the door. Remove all the red coloured parts, they are only used for the transportation of the appliance.
- **When placing, moving, lifting the appliance, do not hold the door handles, do not pull the condenser at the rear part of the refrigerator and do not touch the compressor unit.**
- Remove the two supports **9** from the bag (see Fig. **1**) and insert them into the guides at the top of the back of the refrigerator.
- When installing the refrigerator, do not tilt it more than 30 ° to the intended location.

The appliance **should not be connected** to the mains until all packing and transport materials are removed. Suitably dispose the packaging material.

### First use:

- Wash the inside of the device and all internal parts with lukewarm water and a small amount of ethyl alcohol-based or water-based cleaning agents to remove the new-device smell, and dry well.
- Once the unit has been positioned, wait for about **30 minutes** before it is connected to the mains.


If the transport and installation conditions specified by the manufacturer have not been observed and the appliance has been tilted more than 30° during transport, allow it to stand in an upright position for **at least 4 hours until it is connected to the mains**. Otherwise, the refrigerator compressor may be damaged.

### Storage links:

- Put medicines and other pharmaceutical preparations in the device when the required temperature has been reached inside it.
- Do not cover the circulating air fan inside the cooling slots: Place the drugs or pharmaceuticals in such a way that the air can circulate freely. This is important for unhindered air circulation and uniform temperature distribution throughout the device.
- Refrigerated drugs or pharmaceutical preparations must not touch the back wall.
- Keep liquids in closed containers.

## DESCRIPTION OF THE APPLIANCE, BASIC PARTS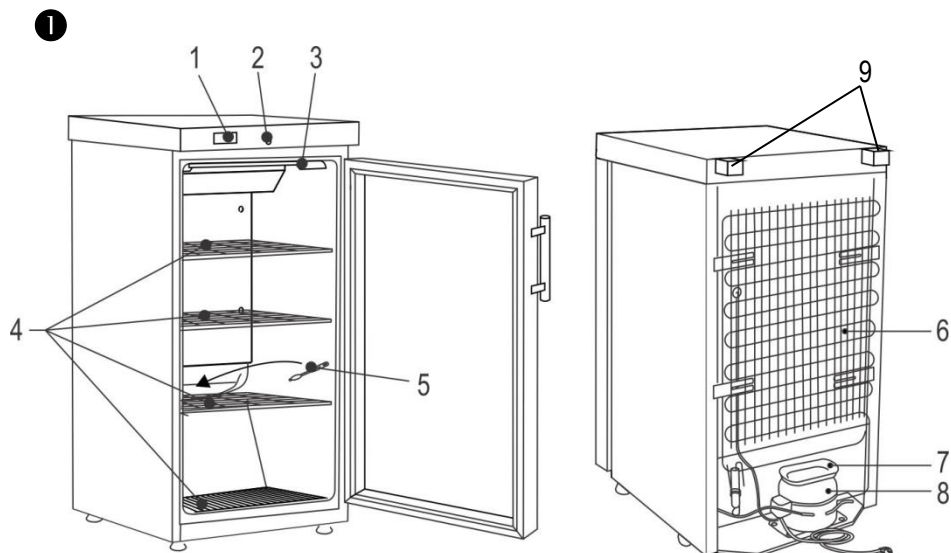

1 Electronic controller

6 Condenser

2 Lock

7 Tray for collection of the defrost water

3 Interior lighting unit

8 Compressor

4 Shelves

9 Distance supports

5 Fan Thaw water channel cleaner

**When the device is turned on, the fan turns on and runs continuously. When the refrigerator door is opened, the fan turns off.**

When the appliance is connected to the mains for the first time, the electronic controller will display the current temperature inside the appliance (hereinafter referred to as the **ACTUAL** temperature).

The average temperature of the device is set to +5 °C and the temperature is not adjustable (The electronic controller display will show the actual temperature range from +2 to +8).

THE SET TEMPERATURE CAN BE CHANGED BY THE MANUFACTURER OR SNAIGĒ SERVISAS OFFICIAL REPRESENTATIVES.

**The electronic controller will always display the actual temperature in the refrigeration unit instead of the set temperature.**

**NOTE.** The electronic controller displays a variable air temperature near the evaporator and is used to control the refrigeration unit. T.y. the actual average temperature is displayed on the controller, so when storing products that are sensitive to temperature changes, it is recommended to use additional temperature measuring devices (arranged at different device heights) for optimal temperature selection.

## ELECTRONIC CONTROLLER

### 1 - On/off button of the refrigeration device. On/off button of lighting

- Press and hold button **1** until **“on”** appears on controller display **3**. Release the button and the refrigerator will start.
- Press and hold button **1** until **“off”** appears on controller display **3**. Release the button and the refrigerator will turn off. **ATTENTION!** The electrical voltage remains unswitched.
- A short press of the button turns the refrigerator lighting on and off.

### 2 - Manual defrost button.

### 3 – Controller display. Meanings of the icons on the screen:

-  lights up on the screen - the compressor is running.
-  lights up on the screen – automatic defrost is in progress.
-  lights up on the screen - the fan inside the device is running.

## Electronic controller indication values

Indication	The reason	What to do?
<b>Hi</b>	The device temperature is too high	Check that the appliance door is closed and tightly closed. If the <b>Hi</b> symbol still flashes after an hour when the door is closed tightly, contact Service.
<b>uHi</b>	Mains voltage higher than 270 V	Mains failure. Contact your electrician.
<b>E01</b>	Temperature sensor (sensor) failure.	Contact Service
<b>dEF</b>	The refrigeration unit is defrosting	Normal operation of the refrigerator. After the defrost time, the refrigerator will operate normally and the display will show the actual temperature.

## CLEANING AND CARE


THE REFRIGERATOR DEFRODS AUTOMATICALLY. The ice droplets formed on the rear wall defrost during the pause of the compressor, at which time the **DEF** indication appears on the electronic controller display. The thawed water drains through the melt water drain to a bath above the compressor (Fig. 1) and evaporates from there.

If, for some reason, the appliance does not defrost automatically, the manual defrost function can be activated: press button **2**, **DEF** will appear on display **3**. The refrigerator compressor will stop freezing: it will run for 20 minutes forced defrosting. After the defrost time, the appliance will start operating normally and the controller display **3** will show the actual temperature inside the appliance.

CLEAN THE REFRIGERATOR REGULARLY.

**Remember that before cleaning, it is mandatory to turn off the refrigerating device with the on/off button and disconnect it from the power supply network - pull the cord plug from the electrical outlet.**

- Protect the appliance inner surfaces and plastic parts from fats, acids and sauces. In case of accidental spills – immediately clean with warm soapy water or dishwashing detergent. Wipe dry.
- Regularly clean the hermetic gaskets of the doors. Wipe dry.
- At least once a year clean dust away from the back part of the appliance body and the compressor. A soft brush, electrostatic cloth or vacuum cleaner may be used for cleaning.
- Regularly clean the thaw water channel space in the refrigerator with a special cleaner for that purpose (see fig.).
- Clean the glass with glass cleaner.
- Clean the inside, shelves and outside of the refrigerator with lukewarm water or ethyl alcohol-based cleaners or water-based household cleaners.
- **DO NOT** use detergents containing abrasive particles, acid, alcohol or benzene for cleaning the appliance inner and outer surfaces. Clean the glass with glass cleaner
- **DO NOT** use for cleaning cloths or sponges that have coarse surfaces intended for scrubbing.
- **DO NOT** use spray ice solvents.
- **DO NOT** clean with compressed air / steam.
- If the appliance is to be left switched off for an extended period switch off the mains voltage, disconnect the battery using the On/Off switch, clean, dry it out and leave its doors open to avoid mould.
- **Only disinfect the device with ethyl alcohol-based products.**

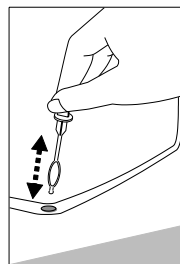## OPERATION PROBLEMS AND THEIR SOLUTIONS

- **The appliance is plugged in to the mains but it does not work.** Check if your house electricity supply installations are in order. Check if the plug is correctly inserted into the electricity supply socket.
- **The noise has become louder.** Check if the appliance is standing stable, on a level place. To make it level, regulate the front feet. Check if the appliance is not touching any furniture and if no part of the refrigeration system at the back part of the appliance body is touching the wall. Pull the appliance away from any furniture or walls. Check if the cause of increased noise isn't due to bottles, cans in the appliance that might be touching each other.

- **Water has appeared at the bottom of the refrigerator compartment.** Check whether a thaw water channel space isn't blocked. Clean the thaw water channel space with a cleaner intended for that purpose.
- **When the appliance door is opened, the rubber sealing gasket pulls out.** The gasket is smeared with sticky foodstuffs (fat, syrup). Clean the sealing gasket and the groove for it with warm water containing soap or dishwashing detergent and wipe dry. Put the rubber sealing gasket back into the place.
- **High temperature in appliance, the compressor operation pauses are short.** Ascertain whether the appliance door/doors are shutting tightly, whether the door wasn't kept open longer than necessary when taking out or putting in foodstuffs or whether a large amount of warm food was placed in the appliance.
- **The top of the appliance has condensation on it.** The ambient relative air humidity is above 75 %. Ventilate the room where the appliance stands and if possible remove the cause of the humidity.
- **The temperature inside the refrigerator exceeds the set permissible limit.** Refrigeration device (evaporator) does not have time to defrost automatically. Activate the manual defrost function: press the button **2**, the indication **dEF** will appear on the display **3**. After the defrost time, the refrigerating appliance will start working in normal mode and the actual temperature inside the appliance will be displayed on the controller display **3**.

**Remarks on the appliance operating noise.** As the appliance operates and goes through the freezing cycle, various noises will be emitted. This is normal and is not a sign of any malfunction. As the refrigerant circulates around the refrigeration system, it causes sounds like murmuring, bubbling, or rustling. Louder sounds like popping or clicking may be heard for a short time as the refrigerator's compressor switches on.

## ENVIRONMENTAL PROTECTION INFORMATION

**Important Note:** This appliance is marked according to the European directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE). By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product.

The symbol on the product, or on the documents accompanying the product, indicates that this appliance may not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment.

For more detailed information about treatment, recovery and recycling of this product, please contact your local authority, your household waste disposal service or the retailer where you purchased the product.

Our packaging are made of environmentally friendly materials, which can be reused:

- The external packaging is made of cardboard/foil
- The FCKW free shape of foamed polystyrene (PS)
- Polyethylene (PE) foils and bags

The refrigerants and foaming agents, which are 100% free of FCKW and FKW have been used for manufacturing our product. Therefore we are helping to protect the ozone layer and prevent the increase of greenhouse effects. Also, the innovative technology and environmentally friendly insulation help in reducing energy consumption.