SNAIGĖ
M E D I C A L

MD35DD-P301CDQ

MD35DD-P600MDQ

MEDICININIS ŠALDYTUVAS LT

Originali naudojimo instrukcija

PHARMACY REFRIGERATOR EN

Original Instruction for use

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DEAR CUSTOMER!

CAREFULLY READ THE INSTRUCTION MANUAL. If instructions are not followed, there is a risk of injury, damage of the device and loss of the right to free warranty service.

The Manufacturer shall not be held liable for damage arising from the failure to observe the instructions contained in this manual.

Keep this manual for future reference, or to pass it over to the next user.

This instruction applies to more than one model and may contain inconsistencies.

General information:

The refrigeration device is intended for the professional storage of medicines, etc. for the storage of pharmaceutical preparations at a temperature of +5 °C degrees.

The device is suitable for medicines, etc. for pharmaceutical preparations that must be stored at a temperature between +2 °C and +8 °C.

The temperature - energy parameters of the device meet the requirements of DIN 13277.

The noise emitted by the device does not exceed 50 dB(A)

Do not use refrigeration inside the device:

- For cooling and storage of chemically unstable, flammable or alkaline substances.
- For cooling and storage of blood, plasma or other infused body fluids.

GUIDELINES CONCERNING THE SAFETY OF USE**General safety requirements:**

- WARNING: Do not cover the ventilation openings of the device.
- WARNING: Do not use electrical devices inside the storage compartments of the device, unless they are of the type recommended by the manufacturer.
- WARNING: Do not use any mechanical tools or other/chemical agents other than those recommended by the manufacturer to speed up the defrosting process.
- WARNING: Do not store explosive substances such as aerosol cans with a flammable propellant in this device.
- WARNING: When positioning the device, ensure the supply cord is not trapped or damaged. This can cause a short circuit and a fire.
- WARNING: Do not locate multiple portable socket-outlets or portable power supplies at the rear of the device.
- WARNING: Disposal of the device should be according to national rules.

WARNING: Do not damage the device refrigeration system. It contains the refrigerant gas R600a. If the refrigeration system is damaged:

- Do not use any open flame.
- Avoid sparks - do not turn on/off any electrical devices or lighting fixtures.
- Immediately ventilate the room: air the room in which the device is placed for a few minutes (the size of the room for a product containing isobutane/ R600a must be at least 4 m³).

C-Pentane is used as blowing agent in the insulation foam and it is a flammable gas.

- This device is designed only for cooling and storing of pharmaceuticals.
- This device is not designed for storage of explosives, such as aerosols with flammable gases.
- It is prohibited to store petrol and other flammable liquids near the device.
- Do not allow children to play with the packaging material from the device.
- Only qualified personnel who are familiar with the installation, commissioning and operation of the refrigerating device are allowed to install, check, maintain and operate the device. Qualified personnel are persons who, due to their technical education, knowledge and experience and knowledge of the relevant standards, are able to assess and perform the work assigned to them and to recognize potential hazards. They must be trained, instructed and authorized to operate the device. Children and persons with limited physical, sensory or mental capabilities or without sufficient experience and knowledge are prohibited from operating and operating this device

Requirements for the connection to the mains:

- The device should ONLY BE CONNECTED TO AN EARTHED SOCKET INSTALLED IN ACCORDANCE WITH THE REGULATIONS. This product complies with all binding CE labelling directives. The earthed electrical socket by which the device is connected to the mains should be in an accessible place. It is a legal requirement that the device is properly earthed. The manufacturer will not be held liable for any damage or injury which may result from the failure to fulfil this requirement.

- It is mandatory to replace a damaged power cord with a proper one of the same type as the one used by the refrigeration device manufacturer to prevent damage to health and property
- To avoid the exposure to danger, always have faulty cable to be replaced only by the manufacturer, by our customer service or by a qualified person and with a cable of the same type.
- It is forbidden to readjust or modify any parts of the unit. It is crucial not to damage the capillary tube visible in the compressor recess. The tube may not be bent, straightened or wound. If the capillary tube is damaged by the user the guarantee will be void.
- Do not use adapters, multiple sockets and two-wire extension leads. If it is necessary to use the extension lead, it must be equipped with a protection ring and a single socket and must have a VDE/GS safety certificate. If an extension lead is used (with a protection ring and safety certificate), its socket must be located at a safe distance, away from the sinks, and must not be in a place where it could be flooded by water or waste water.
- The frequency and power of the electricity supply in your house must conform to the general data parameters of the device as it is shown in the rating label.
- Make sure that the plug of the electrical cord is not squeezed by the back wall of the device or otherwise damaged. **A DAMAGED PLUG CAN BE THE CAUSE OF A FIRE!**

 **CAUTION!** If the installed interior lighting light-emitting diode (LED) lamp for the device is not functioning, contact the maintenance and repair representative for replacements. **LAMPS CAN BE REPLACED ONLY BY THE MAINTENANCE AND REPAIR REPRESENTATIVE**

 **DO NOT TRY REPAIR THE DEVICE ON YOUR OWN. ALWAYS CONTACT AN AUTHORIZED TECHNICIAN.** To avoid the exposure to danger, always have faulty cable be replaced only by the manufacturer, by our customer service or by qualified person with a cable of same type.

Requirements for safe use:

- It is forbidden to use a technically damaged device.
- If the device is unplugged (for cleaning, moving to another place, etc.), it may be repeatedly switched on after **15 min**.
- When performing activities, such as cleaning, maintenance or moving, the device must be completely disconnected from power supply (by pulling the plug out from the socket). (Do not forget to turn off the emergency (reserve) power supply of the electronic controller with button **11** (see fig.1) before that.
- **Do not pull on the cord, but hold in the body of the plug instead.**
- Do not place heavy objects on the top of the refrigerating device.
- Do not place any switched on electrical devices on top of the device because this may cause ignition of plastic parts.
- Do not place any dishes with liquids on top of the device.
- Do not climb on or sit on the device, do not lean on or hang on the device doors.
- Do not use water spray and steam to clean the device.

WARNING: the device may only be used with original accessories from the manufacturer or accessories from other suppliers approved by the manufacturer. Use of non-approved accessories is at the user's own risk.

It is recommended to transport the refrigerating device in a vertical position; THE MANUFACTURER IS NOT RESPONSIBLE FOR FAILURES OF THE DEVICE IF THE CONDITIONS OF TRANSPORTATION NOT IS FOLLOWED.

INSTALLATION AND OPERATING CONDITIONS


Positioning:

- This device can be installed in a dry, well ventilated indoor location where the ambient temperature corresponds to the climate class indicated on the rating plate of the device.

CAUTION! Do not install this device in areas that are too humid or too cold. Place the device away from heat sources such as radiators, or direct sunlight. The device must not touch any pipes for heating, gas or water supply, or any other electrical devices.

- Do not cover the ventilation holes at the top of the devices – it must be a good air circulation around the device. There should be a gap of at least 10 cm between the top of the device body and any furniture that may be above it. If this requirement is not followed, the device consumes more electrical energy and its compressor may overheat.
- The device must stand on a level surface and must not touch the wall. If necessary, regulate the height of the device by adjusting the levelling feet: by turning them clockwise – the front of the device rises, by turning them counter clockwise – it comes down. If the device is tilted slightly backward – the doors will close by themselves.
- **Scheme of the external space required to operate the device. External dimensions for installation.** If the device is placed in a corner, a gap of at least 5 cm must be left between the device body and the wall (see Fig.).

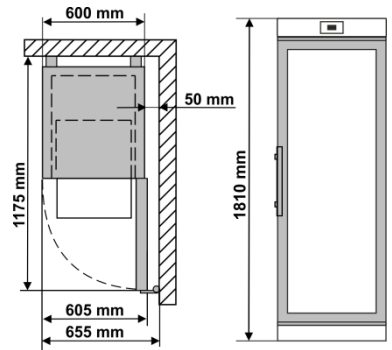
Preparing for operation:

- It is recommended to prepare the device for operation with a helper.
- Remove package. Lift the device away from foamed polystyrene base. Therefore take out adhesive tapes from sides and simply lift the device up and take out the board. Strip adhesive tapes off and open the door. Remove all the red coloured parts, they are only used for the transportation of the device.
- When positioning the device in chosen location, it will move more easily into position if you lift the front a little and incline it backward allowing it to roll on its casters.

When placing, moving, lifting the device, do not hold the door handles, do not pull the condenser at the rear part of the refrigerator and do not touch the compressor unit. When placing the device in its intended location, it is not recommended to turn it over

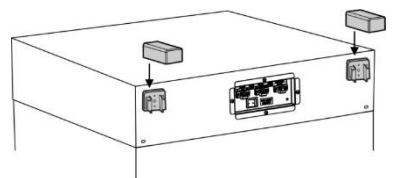- Remove the handle, the screws and tools from the plastic bag. Screw the handle with screws in the holes prepared on the side of the door.
- Remove the two supports from the bag and insert them into the guides at the top of the back of the device (see fig.)

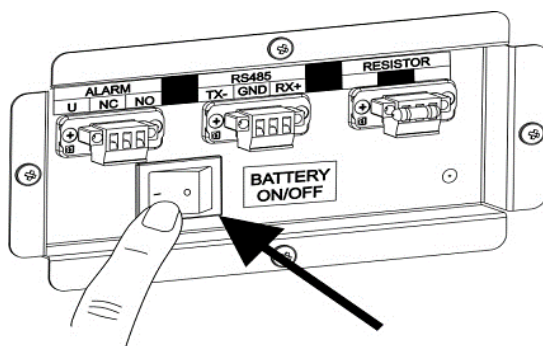
 The device **should not be connected** to the mains until all packing and transport materials are removed. Suitably dispose the packaging material.

First use:

- Wash the interior of the device and all internal parts with lukewarm water and a small amount of ethyl alcohol-based cleaning agents - this will remove the characteristic smell of a new device - and dry well
- Once the unit has been positioned, wait for about **30 minutes** before it is connected to the mains.

 After connecting the device to the electrical network, use button **11** (see fig. 1) to turn on the emergency (reserve) power supply of the electronic controller. A symbol  appears on the controller screen.

 When the device is decommissioned and disconnected from the mains, it is necessary to disconnect the emergency backup power supply to the controller using switch **11** (Fig. 1)

 **If the unit has been tilted more than 30° during transport: Wait at least 4 hours before connecting the device to the power supply. If this is not done, the device compressor may fail.**

Storage links:


- Put medicines and other pharmaceutical preparations in the device when the required temperature has been reached inside it.

 Do not cover the circulating air fan inside the cooling slots. **Place the drugs or pharmaceuticals in such a way that the air can circulate freely.** This is important for unhindered air circulation and uniform temperature distribution throughout the device.

- Refrigerated drugs or pharmaceutical preparations must not touch the back wall.
- Keep liquids in closed containers.

DESCRIPTION OF THE DEVICE, BASIC PARTS


Fig. 1

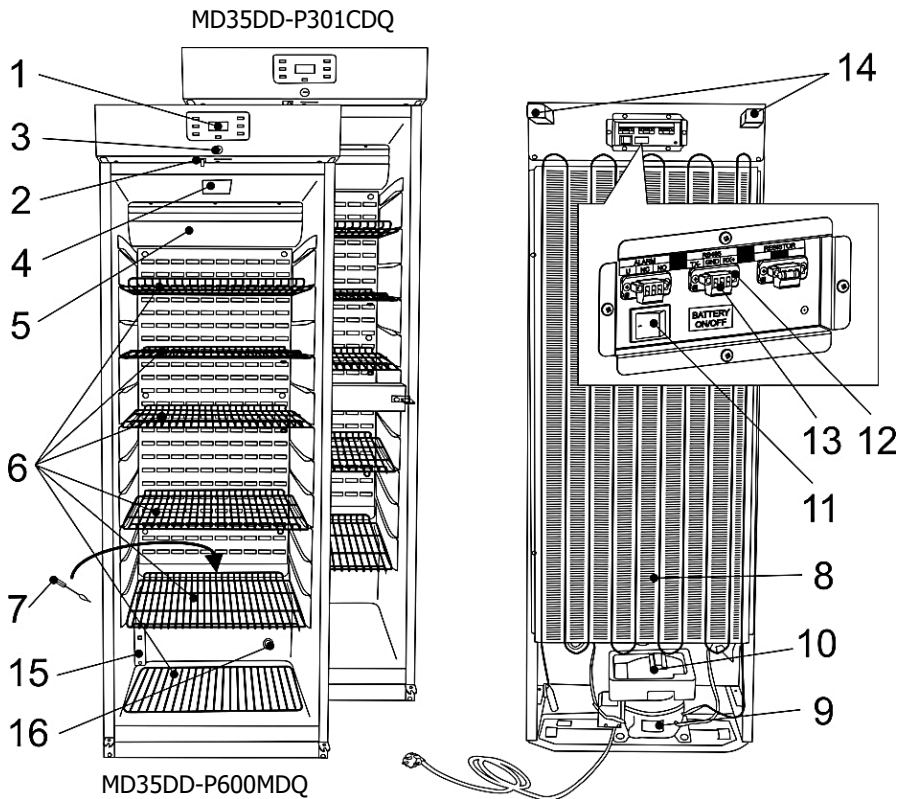1	Electronic controller	8	Condenser
2	Fan and interior lighting switch	9	Compressor
3	Lock*	10	Tray for collection of the defrost water
4	Interior lighting	11	Emergency power disconnect switch
5	Fan	12	External alarm connection
6	Shelves	13	Network connection RS 485
7	Thaw water channel cleaner		

* Lock keys (2 pcs.) are attached to one of the shelves.

Maximum load per each shelf is 40 kg

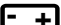After connecting the device to the mains, switch on the emergency backup power supply to the controller using switch **11** (Fig. 1) A symbol  appears on the controller screen **When the device is turned on, the fan turns on and runs continuously. When the device door is opened, the fan turns off and the interior lighting of the device turns on.**

The lighting is controlled by an electronic controller. If needed, the interior lighting can be turned on for continuous operation, press the button 

THE DEVICE HAS A SECOND SAFETY CONTROL THAT PREVENTS THE TEMPERATURE INSIDE THE DEVICE FROM FALLING BELOW +2 °C.

When the device is connected to the mains for the first time, the electronic controller will display the current temperature inside the appliance (hereinafter referred to as the **ACTUAL** temperature).

 **The electronic controller will always display the actual temperature in the device instead of the set temperature.**

The average temperature of the device is set to +5 °C and the temperature is not adjustable (+5 °C to ensure an average temperature, the electronic controller display will show the actual temperature range from +2 to +8). THE SET TEMPERATURE CAN BE CHANGED BY THE MANUFACTURER OR OFFICIAL REPRESENTATIVES.

NOTE. The electronic controller on the display shows the variable air temperature near the evaporator and is used to control the refrigeration unit. I.e. the controller displays the actual average temperature, so **when storing medications that are sensitive to temperature changes**, it is recommended to use additional temperature measuring devices (positioned at different device heights) for optimal temperature selection.

Technical data sheet

Model	MD35DD-P301CDQ	MD35DD-P600MDQ
Gross capacity, l	350	
Nett capacity, l	230	
External dimensions (H x W x D),mm	1810 x 600 x 620	
Weight: Brutto / netto, kg	82 / 77	82 / 77
Rated current, A	1,7	
Voltage / Frequency rated, V / Hz	220 – 240 / 50	
Rated power input, W	230	
Climate class	7 (nuo +10 iki +35, ≤75%)	
Product temperature range (Min/Max), °C	nuo +2 iki +8	
Annual energy consumption, kWh/year	648	648

DESCRIPTION OF THE ELECTRONIC CONTROLLER, TEMPERATURE ADJUSTMENT

	Pressing the SET button displays the set temperature. In programming mode it selects a parameter or confirm an operation
	Pressing the UP button takes you to the quick menu. In programming mode, scroll through parameter codes or increase the displayed value.
	Pressing the DOWN button in programming mode it browses the parameter codes or decreases the displayed value.
	DATA exporting data to a USB storage device
	Pressing the button turns on the interior lighting.
	REC - saving data in the memory of the electronic controller. The data save function is activated at the factory. To disable this function, you need to contact the service department.

Electronic controller display symbols and their explanation

Symbol	Operation	Rated
	On	Compressor enabled
	Flashing	Compressor operation protection activated.
	On	The automatic defrost function is activated
	Flashing	Steam drying is in progress
	On	Compressor enabled
	Flashing	Po automatinio atitirpinimo trumpam sustabdytas ventiliatoriaus veikimas. Normalus šaldymo prietaiso veikimas.
	On	Critical errors are logged.
	On	Temperature recording is activated.
	On	Battery works OK.
	Flashing	BATTERY NOT CHARGING: The emergency power supply is not switched on (see Fig. 1, switch 11) or a fault due to battery charging / failure.
°C/°F	On	Measurement unit
	Flashing	Programming phase
	On	The interior lighting enabled

Set the date and time settings, disable the real-time menu (the RTC) alarm.

<ol style="list-style-type: none"> 1. Enter the Programming Mode pressing both SET+DOWN keys for 3 sec. 2. Use the UP or DOWN button till the CLO menu is displayed
Press SET to enter the settings menu
<p>HUr = hour → press SET to modify/confirm. UP/DOWN to modify the parameter.</p> <p>Min = minutes → press SET to modify/confirm. UP/DOWN – to modify the parameter.</p> <p>dAy = Day of the month → press SET to modify/confirm. UP/DOWN – to modify the parameter.</p> <p>Udy = Day of the week → (Sun, Mon, Tue, Ued,thE, Fri, Sat) press SET to modify/confirm. UP/DOWN – to modify the parameter.</p> <p>Mon = month → press SET to modify/confirm. UP/DOWN – to modify the parameter.</p> <p>yEA =year → press SET to modify/confirm. UP/DOWN – to modify the parameter.</p>
Press both SET and UP together The audible alarm is switched off at the same time. After not pressing any button for about 10 s. automatically exits RTC.

 **When the refrigerator is working in emergency mode (in case of power failure), only the temperature change inside the device is registered. Compressor, fan, lighting do not work.**

Temperature recording and data transfer to a USB key

To transfer data from the electronic controller to a USB stick, follow these steps:

1. Insert the supplied product-specific USB storage media that comes with the device;
2. Press the **DATA** button until **dat** appears;
3. The code **Lod** will be displayed on the screen during data export, and then a code **Alr**;
4. When the data transfer is complete, the following code will be displayed:
 - **End** if everything's okay
 - **Err** if the data transfer failed.

The last ones are stored in the memory of the electronic controller:

- a) 100 alarm entries along with the date and duration of the event „ALARM.txt”
- b) A summary of temperature measurement records along with the date and time of the recording „DATA.txt.”

CLEANING AND CARE


THE DEVICE DEFROSTS AUTOMATICALLY. The ice droplets formed on the wall of the evaporator are defrosted during the pause of the compressor operation, at which time the automatic defrost sign ❄️ is activated on the display of the electronic controller. The thawed water drains through a tube into the bath **10** (Fig. 1), designed to collect the melt water, and evaporates from there.

CLEAN THE DEVICE REGULARLY.

Remember that before cleaning, it is necessary to turn off the emergency (reserve) power supply of the electronic controller with button 11 (fig. 1), and to disconnect it from the power supply network - pull the cord plug from the electrical outlet.

- Regularly clean the hermetic gaskets of the doors. Wipe dry.
- REGULARLY CLEAN THE THAW WATER CHANNEL SPACE IN THE device WITH A SPECIAL CLEANER FOR THAT PURPOSE (see fig.).
- At least once a year, clean the dust accumulated on the back part of the refrigerating device body, the condenser and the compressor. You can use a soft brush, electrostatic wipes or a vacuum cleaner for cleaning.
- Clean the inside, shelves and outside of the device with lukewarm water or ethyl alcohol-based cleaners or water-based household cleaners.
- Clean the glass with products intended for glass.
- DO NOT use cleaners containing abrasive particles, acid, gasoline to clean the inside and outside of the device, the door seal gasket.
- DO NOT use cloths or sponges for cleaning on rough, scrubbing surfaces.
- DO NOT use spray deicers.
- DO NOT clean with compressed air/steam.
- If the device is left empty for a long time, turn it off, defrost, clean, dry and leave the door open to prevent mold from forming inside the appliance.
- **Only disinfect the device with ethyl alcohol-based products.**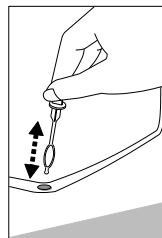

CHANGING THE OPENING DIRECTION OF THE DOORS

We recommend changing the opening direction of the device door with an assistant. You will need: wrench 8mm and 10mm and a Phillips screwdriver.

When changing the door opening direction, it is FORBIDDEN to place the device horizontally.

Before you start, please note the following:

- Always ensure that the correct number of washers is used on all hinge bolts and screws.
- So that the doors can close tightly against the frame, take the utmost care that they are properly aligned during reassembly.

ATTENTION! Switch off the device and remove the plug from the power supply socket. Do not forget to turn off the emergency (reserve) power supply of the electronic controller beforehand.

For MD35DD-P301CDQ

Do the actions below in the order they are listed:

1. Turn the device at a 45° angle to the back and support it firmly

2. Unscrew the two screws **7** and remove the lower bracket **8** together with the set of washers **6** on the axle and the spacer **5**.

3. Unscrew the two screws and remove the bracket **10**.

4. Remove the bottom door **18**.

5. Unscrew the fixing screws **17** of the middle bracket and remove the middle bracket **16** together with the washers on the axle and the spacer **15**.

6. Remove the upper door **4**.

7. Unscrew the axle **2** together with the nut **1** and the washers **3** from the upper bracket and screw the whole set on the opposite side of the refrigerator;

8. Remove the blinds **19** in the cross section of the cabinet and install them on the opposite side of the cabinet.

9. Unscrew the fastening screws **12** from both doors, remove the handles **11**. Pull out the blinds **13** and **14** and install them on the opposite side of the door. Fix handles **11** in the loosened holes with screws **12**.


10. Hook the upper door **4** after inserting the upper hinge into the door.

11. Screw the middle bracket **16** with a set of washers and gaskets on the opposite side of the cabinet, before inserting the bracket's shoulder into the upper door.

12. Hook the bottom door **18** on the axle of the middle bracket.

13. Plate **9** remains in place. Screw the bottom bracket **8** with a set of washers and spacers on the opposite side of the cabinet, before inserting the bracket's shoulder into the door.

14. Screw the bracket **10** on the opposite side of the cabinet.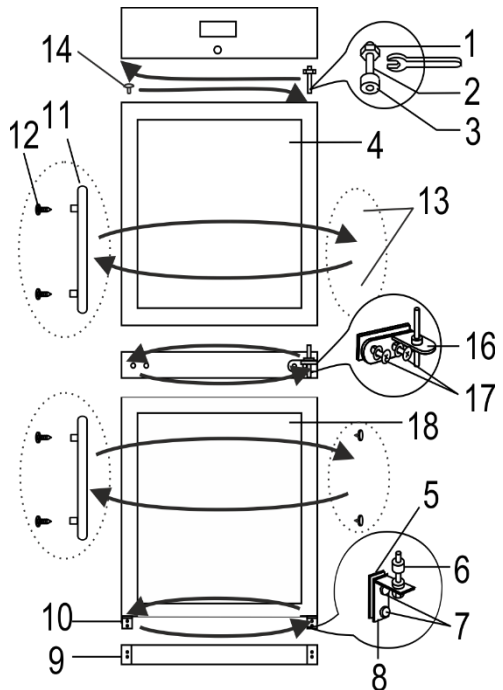

For MD35DD-P300MDQ

Do the actions below in the order they are listed:

1. Turn the device at a 45° angle to the back and support it firmly

2. Unscrew the two screws **7** and remove the lower bracket **8** together with the set of washers **6** on the axle and the spacer **5**.

3. Unscrew the two screws **7** and remove the bracket **10**.

4. Remove door **4**.

5. Unscrew the axle **2** together with the nut **1** and the washers **3** from the upper bracket and screw the whole set on the opposite side of the refrigerator.

6. Hang the door after turning it 180° and inserting the upper hinge into the door.

7. Remove the blind **13** from the door, unscrew the fixing screws **12**, remove the handle **11** and fix it by screwing the screws into the free holes on the same side. Insert the blind **13** into the remaining hole.


8. Plate **9** remains in place. Screw the bottom bracket **8** with a set of washers and spacers on the opposite side of the cabinet, before inserting the bracket's shoulder into the door.

9. Screw the bracket **10** on the opposite side of the cabinet.

If necessary, adjust the even position and fit of the door:

- adjust the even position of the door by sliding the lower bracket left / right;
- Ensure that the door fits snugly by changing the amount of spacers under the brackets.

NOTE. If you tilt the device at an angle of more than 45 ° when changing the direction of the door opening, allow it to stand upright for 4 hours before connecting it to the mains.

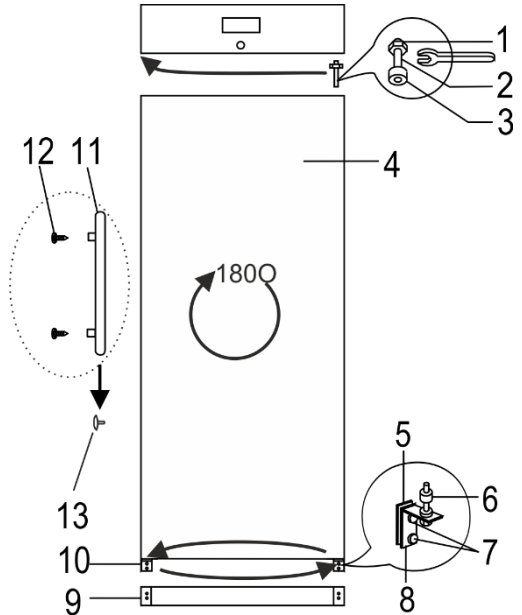
HOW TO USE THE ACCESS PORT

You will need 4mm flat head screwdriver.

ATTENTION! Switch off the device and remove the plug from the power supply socket. Do not forget to turn off the emergency (reserve) power supply of the electronic controller beforehand.

1. Remove the sealing paste from the access port hole on side **A** of the compressor compartment (see Fig. 1).
2. Inside the device **B**, remove the cover from the access port hole using a screwdriver (see Fig. 2).
3. Push out the insulating material with a screwdriver (see Fig. 3).
4. Insert the sensor cable from inside the device B through the hole (see Fig. 4).
5. On the niche side A, wrap the sensor wire with insulating material and push it into the hole of the access port (see Fig. 5).
6. Additionally, seal the hole on side A with sealing paste (see Fig. 6).

After connecting the device to the electrical network, to turn on the emergency (reserve) power supply of the electronic controller.

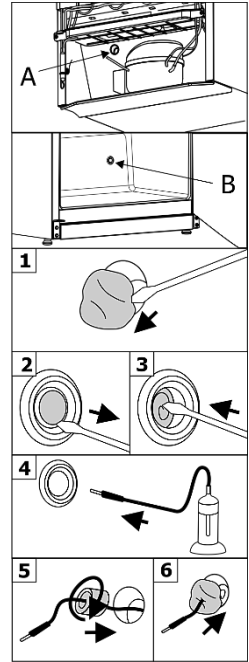OPERATION PROBLEMS AND THEIR SOLUTIONS

- **The appliance is plugged in to the mains but it does not work.** Check if your house electricity supply installations are in order. Check if the plug is correctly inserted into the electricity supply socket.
- **The symbol  flashes. The battery is not charging.** Check that the emergency power supply is switched on (detailed description in the INSTALLATION AND PREPARATION section). If the display does not go out when the emergency power is switched on, the fault is due to the battery being charged / faulty. Contact Service.
- **The noise has become louder.** Check if the appliance is standing stable, on a level place. To make it level, regulate the front feet. Check if the appliance is not touching any furniture and if no part of the refrigeration system at the back part of the appliance body is touching the wall. Pull the appliance away from any furniture or walls. Check if the cause of increased noise isn't due to bottles, cans in the appliance that might be touching each other.
- **Water has appeared at the bottom of the refrigerator compartment.** Check whether a thaw water channel space isn't blocked. Clean the thaw water channel space with a cleaner intended for that purpose.
- **When the appliance door is opened, the rubber sealing gasket pulls out.** The gasket is smeared with sticky foodstuffs (fat, syrup). Clean the sealing gasket and the groove for it with warm water containing soap or dishwashing detergent and wipe dry. Put the rubber sealing gasket back into the place.
- **The top of the appliance has condensation on it.** The ambient relative air humidity is above 75 %. Ventilate the room where the appliance stands and if possible remove the cause of the humidity.

- **Alarms codes.** The alarm code is displayed as long as the alarm condition persists. All alarm codes are displayed alternately with the internal temperature, except for **PF1**, which flashes. The visual indication of all faults is accompanied by an audible signal. This audio signal can be turned off by pressing any button. Device has the possibility to connect an external emergency alarm.

Code	The reason	What to do?
PF1	Defective control sensor	Contact Service
PF2	Defective evaporator sensor	Contact Service
PF3	Defective temperature registration sensor	Contact Service
HA1	High temperature	Check operation of the device, ambient temperature.
HA3	High product temperature	Check operation of the device, ambient temperature.
LA1	Low temperature	Contact Service
dA	Open door alarm	Close the door tightly
CA	Severe alarm	Contact Service
EA	External alarm	Contact Service

Notes on the operation sound of the refrigerating device. When the device is working, various operating sounds are emitted during the refrigerating process, it is a normal phenomenon, it is not a malfunction. Hissing, bubbling, hissing sounds are caused by the refrigerant circulating in the refrigeration system. The compressor of the device emits a stronger humming or clicking sound when it starts up for a short time.

ENVIRONMENTAL PROTECTION INFORMATION

Important Note:

This device is marked according to the European directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE). By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product.

The symbol on the product, or on the documents accompanying the product, indicates that this device may not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment.

■ A rechargeable battery is installed in the device. The battery must be removed when you wish to dispose of the device and sent for separate waste treatment for batteries. **Do not damage or short circuit the battery.**

Disposal must be carried out in accordance with local environmental regulations for waste disposal. For more detailed information about treatment, recovery and recycling of this product, please contact your local authority, your household waste disposal service or the retailer where you purchased the product.

Our packaging are made of environmentally friendly materials, which can be reused:

- The external packaging is made of cardboard/foil
- The FCKW free shape of foamed polystyrene (PS)
- Polyethylene (PE) foils and bags

The refrigerants and foaming agents, which are 100% free of FCKW and FKW have been used for manufacturing our product. Therefore we are helping to protect the ozone layer and prevent the increase of greenhouse effects. Also, the innovative technology and environmentally friendly insulation help in reducing energy consumption.

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AB SNAIGĖ