**SNAIGĖ**  
M E D I C A L

**MD40DC-P301BEQ**

**MD40DC-P600BEQ**

**MEDICININIS ŠALDYTUVAS LT**

Originali naudojimo instrukcija

**PHARMACY REFRIGERATOR EN**

Original Instruction for Use

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**DEAR CUSTOMER!**

CAREFULLY READ THE INSTRUCTION MANUAL. If instructions are not followed, there is a risk of injury, damage of the device and loss of the right to free warranty service.

The Manufacturer shall not be held liable for damage arising from the failure to observe the instructions contained in this manual.

Keep this manual for future reference, or to pass it over to the next user.

This instruction applies to more than one model and may contain inconsistencies.

**General information:**

The refrigeration device is intended for the professional storage of medicines, etc. for the storage of pharmaceutical preparations at a temperature of +5 °C degrees.

The device is suitable for medicines, etc. for pharmaceutical preparations that must be stored at a temperature between +2 °C and +8 °C.

The temperature - energy parameters of the device meet the requirements of DIN 13277.

The noise emitted by the device does not exceed 44 dB(A)

**Do not use refrigeration inside the device:**

- For cooling and storage of chemically unstable, flammable or alkaline substances.
- For cooling and storage of blood, plasma or other infused body fluids.

**GUIDELINES CONCERNING THE SAFETY OF USE****General safety requirements:**

- WARNING: Do not cover the ventilation openings of the device.
- WARNING: Do not use electrical devices inside the storage compartments of the device, unless they are of the type recommended by the manufacturer.
- WARNING: Do not use any mechanical tools or other/chemical agents other than those recommended by the manufacturer to speed up the defrosting process.
- WARNING: Do not store explosive substances such as aerosol cans with a flammable propellant in this device.
- WARNING: When positioning the device, ensure the supply cord is not trapped or damaged. This can cause a short circuit and a fire.
- WARNING: Do not locate multiple portable socket-outlets or portable power supplies at the rear of the device.
- WARNING: Disposal of the device should be according to national rules.

**WARNING:** Do not damage the device refrigeration system. It contains the refrigerant gas R600a. If the refrigeration system is damaged:

- Do not use any open flame.
- Avoid sparks - do not turn on/off any electrical devices or lighting fixtures.
- Immediately ventilate the room: air the room in which the device is placed for a few minutes (the size of the room for a product containing isobutane/ R600a must be at least 4 m<sup>3</sup>).

C-Pentane is used as blowing agent in the insulation foam and it is a flammable gas.

- This device is designed only for cooling and storing of pharmaceuticals.
- It is prohibited to store petrol and other flammable liquids near the device.
- Do not allow children to play with the packaging material from the device.
- Only qualified personnel who are familiar with the installation, commissioning and operation of the refrigerating device are allowed to install, check, maintain and operate the device. Qualified personnel are persons who, due to their technical education, knowledge and experience and knowledge of the relevant standards, are able to assess and perform the work assigned to them and to recognize potential hazards. They must be trained, instructed and authorized to operate the device. Children and persons with limited physical, sensory or mental capabilities or without sufficient experience and knowledge are prohibited from operating and operating this device

### **Requirements for the connection to the mains:**

- The device should **ONLY BE CONNECTED TO AN EARTHED SOCKET INSTALLED IN ACCORDANCE WITH THE REGULATIONS**. This product complies with all binding CE labelling directives. The earthed electrical socket by which the device is connected to the mains should be in an accessible place. It is a legal requirement that the device is properly earthed. The manufacturer will not be held liable for any damage or injury which may result from the failure to fulfil this requirement.
- It is mandatory to replace a damaged power cord with a proper one of the same type as the one used by the refrigeration device manufacturer to prevent damage to health and property
- To avoid the exposure to danger, always have faulty cable to be replaced only by the manufacturer, by our customer service or by a qualified person and with a cable of the same type.

- It is forbidden to readjust or modify any parts of the unit. It is crucial not to damage the capillary tube visible in the compressor recess. The tube may not be bent, straightened or wound. If the capillary tube is damaged by the user the guarantee will be void.
- Do not use adapters, multiple sockets and two-wire extension leads. If it is necessary to use the extension lead, it must be equipped with a protection ring and a single socket and must have a VDE/GS safety certificate. If an extension lead is used (with a protection ring and safety certificate), its socket must be located at a safe distance, away from the sinks, and must not be in a place where it could be flooded by water or waste water.
- The frequency and power of the electricity supply in your house must conform to the general data parameters of the device as it is shown in the rating label.
- When positioning the device, be careful that the electrical cord isn't squeezed in order to avoid damaging it. Do not store heavy objects such as cooling devices or other devices next the device in such a way that they could squeeze and damage the electrical cord. This can cause a short circuit and a fire.
- Make sure that the plug of the electrical cord is not squeezed by the back wall of the device or otherwise damaged. **A DAMAGED PLUG CAN BE THE CAUSE OF A FIRE!**

**CAUTION!** If the installed interior lighting light-emitting diode (LED) lamp for the cooling device is not functioning, contact the maintenance and repair representative for replacements. **LAMPS CAN BE REPLACED ONLY BY THE MAINTENANCE AND REPAIR REPRESENTATIVE**

**DO NOT TRY REPAIR THE DEVICE ON YOUR OWN. ALWAYS CONTACT AN AUTHORIZED TECHNICIAN. TO AVOID THE EXPOSURE TO DANGER, ALWAYS HAVE FAULTY CABLE BE REPLACED ONLY BY THE MANUFACTURER, BY OUR CUSTOMER SERVICE OR BY QUALIFIED PERSON WITH A CABLE OF SAME TYPE.**

### **Requirements for safe use:**

- It is forbidden to use a technically damaged device.
- If the device is unplugged (for cleaning, moving to another place, etc.), it may be repeatedly switched on after **15 min**.
- When performing activities, such as cleaning, maintenance or moving, the device must be completely disconnected from power supply (by pulling the plug out from the socket).
- **Do not pull on the cord, but hold in the body of the plug instead.**
- Do not place heavy objects on the top of the refrigerating device.

- Do not place any switched on electrical devices on top of the device because this may cause ignition of plastic parts.
- Do not place any dishes with liquids on top of the device.
- Do not climb on or sit on the device, do not lean on or hang on the device doors.
- Do not use water spray and steam to clean the device.

**WARNING:** the device may only be used with original accessories from the manufacturer or accessories from other suppliers approved by the manufacturer. Use of non-approved accessories is at the user's own risk.

 **It is recommended to transport the refrigerating device in a vertical position;** THE MANUFACTURER IS NOT RESPONSIBLE FOR FAILURES OF THE DEVICE IF THE CONDITIONS OF TRANSPORTATION NOT IS FOLLOWED.

## INSTALLATION AND OPERATING CONDITIONS

### Positioning:


- This device can be installed in a dry, well ventilated indoor location where the ambient temperature corresponds to the climate class indicated on the rating plate of device.

 **CAUTION!** Do not install this device in areas that are too humid or too cold. Place the device away from heat sources such as radiators, or direct sunlight. The device must not touch any pipes for heating, gas or water supply, or any other electrical devices.

- Do not cover the ventilation holes at the top of the devices – it must be a good air circulation around the device. There should be a gap of at least 10 cm between the top of the device body and any furniture that may be above it. If this requirement is not followed, the device consumes more electrical energy and its compressor may overheat.


- The device must stand on a level surface and must not touch the wall. If necessary, regulate the height of the device by adjusting the levelling feet: by turning them clockwise – the front of the device rises, by turning them counter clockwise – it comes down.

- **Scheme of the external space required to operate the device. External dimensions for installation.** If the device is placed in a corner, a gap of at least 5 cm must be left between the device body and the wall (see Fig.).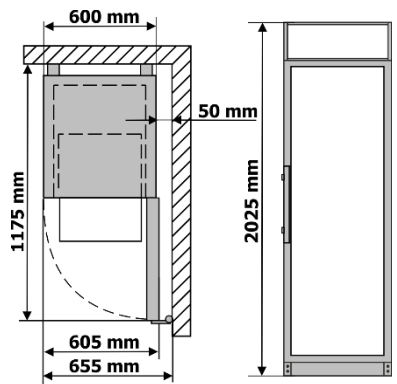

## Preparing for operation:

- It is recommended to prepare the device for operation with a helper.
- Remove package. Lift the device away from foamed polystyrene base. Therefore take out adhesive tapes from sides and simply lift the device up and take out the board. Strip adhesive tapes off and open the door. Remove all the red coloured parts, they are only used for the transportation of the device.
- When positioning the device in chosen location, it will move more easily into position if you lift the front a little and incline it backward allowing it to roll on its casters.
- **When placing, moving, lifting the device, do not hold the door handles, do not pull the condenser at the rear part of the refrigerator and do not touch the compressor unit.** When placing the device in its intended location, it is not recommended to turn it over.
- Remove the handle, the screws and tools from the plastic bag. Screw the handle with screws in the holes prepared on the side of the door.
- Remove the two supports from the bag and insert them into the guides at the top of the back of the device (see fig.)

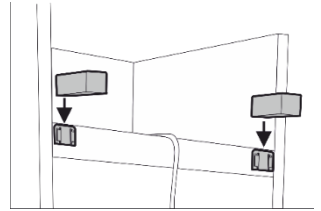 The device **should not be connected** to the mains until all packing and transport materials are removed. Suitably dispose the packaging material.

## First use:

- Wash the interior of the device and all internal parts with lukewarm water and a small amount of ethyl alcohol-based cleaning agents - this will remove the characteristic smell of a new device - and dry well
- Once the unit has been positioned, wait for about **30 minutes** before it is connected to the mains.

After connecting the device to the mains, press and hold **On/off** button until **on** appears on screen. Release the button and the device will start.


 **If the unit has been tilted more than 45° during transport:** Wait at least 4 hours before connecting the device to the power supply. If this is not done, the device compressor may fail.

## Storage links:

- Put medicines and other pharmaceutical preparations in the device when the required temperature has been reached inside it.
-  Do not cover the circulating air fan inside the cooling slots. **Place the drugs or pharmaceuticals in such a way that the air can circulate freely.** This is important for unhindered air circulation and uniform temperature distribution throughout the device.
- Refrigerated drugs or pharmaceutical preparations must not touch the back wall.
- Keep liquids in closed containers.

# DESCRIPTION OF THE DEVICE, BASIC PARTS

**Fig. 1**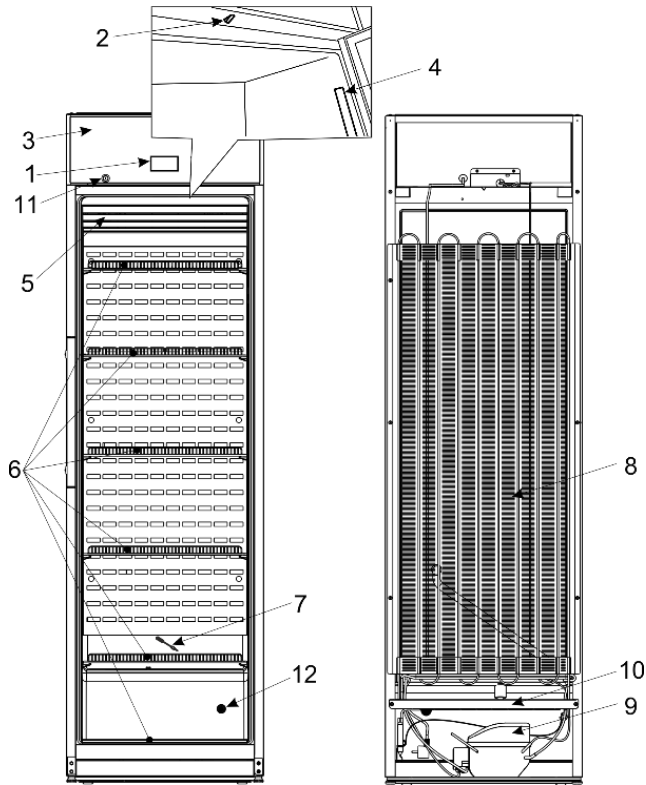

<b>1</b> Electronic controller	<b>7</b> Thaw water channel cleaner
<b>2</b> Fan and interior lighting switch	<b>8</b> Condenser
<b>3</b> Panel	<b>9</b> Compressor
<b>4</b> Interior lighting	<b>10</b> Tray for collection of the defrost water
<b>5</b> Fan	<b>11</b> Lock*
<b>6</b> Shelves	<b>12</b> Access port

\* Lock keys (2 pcs.) are attached to one of the shelves.

Maximum load per each shelf is 40 kg.

**When the device is turned on, the fan turns on and runs continuously. When the device door is opened, the fan turns off and the interior lighting of the device turns on.** If needed, the interior lighting can be turned on for continuous operation, a short press of the **On/off** button.

When the device is connected to the mains for the first time, the electronic controller will display the current temperature inside the device (hereinafter referred to as the **ACTUAL** temperature).

 **The electronic controller will always display the actual temperature in the refrigeration unit instead of the set temperature.**

The average temperature of the device is set to +5 °C and the temperature is not adjustable (+5 °C to ensure an average temperature, the electronic controller display will show the actual temperature range from +2 to +8 °C).


A fixed unregulated control temperature of +5 °C is programmed in the controller. THE SET TEMPERATURE CAN BE CHANGED BY THE MANUFACTURER OR OFFICIAL REPRESENTATIVES.

**NOTE.** The electronic controller displays a variable air temperature near the evaporator and is used to control the device. T.y. the actual average temperature is displayed on the controller, so when storing products that are sensitive to temperature changes, it is recommended to use additional temperature measuring devices (arranged at different device heights) for optimal temperature selection.

## Technical data sheet

Model	MD40DC-P301BEQ	MD40DC-P600BEQ
Gross capacity, l	385	
Nett capacity, l	290	
External dimensions (H x W x D),mm	2025 x 600 x 620	
Weight: Brutto / netto, kg	82 / 76,5	
Rated current, A	1,5	
Voltage / Frequency rated, V / Hz	220 – 240 / 50	
Rated power input, W	250	
Climate class	7 (from +10 to +35 °C , ≤75%)	
Product temperature range (Min/Max), °C	from +2 to +8	
Annual energy consumption, kWh/year	467	348

# ELECTRONIC CONTROLLER

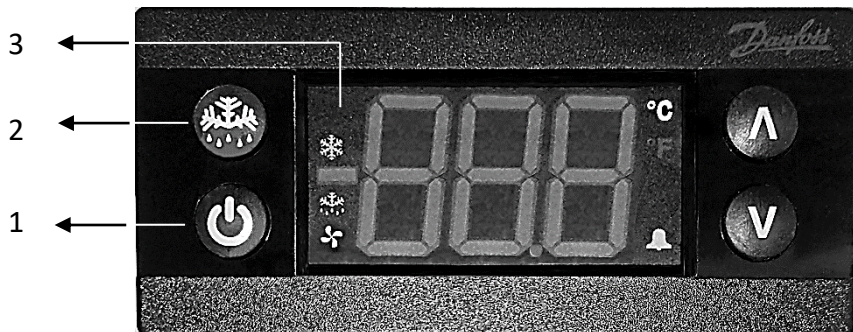## 1 - On/off button of the device.

- Press and hold button **1** until **on** appears on screen **3**. Release the button and the device will start.
- Press and hold button **1** until **off** appears on screen **3**. Release the button and the device will turn off. **ATTENTION!** The electrical voltage remains unswitched.
- A short press of the button turns the device lighting on and off.

**NOTE.** In a device with an solid door does not control the lighting. The lighting turns on automatically when the door is opened and turns off when the door is closed.

## 2 - Manual defrost button.

**3 – Controller display.** Meanings of the icons on the screen:

	lights up on the screen - the compressor is running.
	lights up on the screen – automatic defrost is in progress.
	lights up on the screen - the fan inside the device is running.
	lights up on the screen – emergency temperature change or device failure. When the alarm is active, a 10-second beep sounds periodically every minute. The alarm will turn off when normal working conditions are reached in the device.
“ . ”	Only the segment lights up on the screen - the controller is disabled.

## Electronic controller indication values

Indication	The reason	What to do?
<b>Hi</b>	The temperature in the device is too high	Check that the device door is closed and tightly closed. If the <b>Hi</b> symbol still flashes after an hour when the door is closed tightly, contact Service.
<b>Lo</b>	The temperature in the device is too low	Contact Service
<b>uHi</b>	Mains voltage higher than 270 V	Mains failure. Contact your electrician.
<b>E01</b>	Temperature sensor (sensor) failure.	Contact Service
<b>E02</b>	Temperature sensor (sensor) failure.	Contact Service
<b>dor</b>	The door is not closed	Close the door.
<b>def</b>	The device is defrosting	Normal operation of the device. After the defrost time, the refrigerator will operate normally and the display will show the actual temperature.

## CLEANING AND CARE


THE DEVICE DEFRODS AUTOMATICALLY. The ice droplets formed on the rear wall defrost during the pause of the compressor, at which time the **def** indication appears on the electronic controller display. The thawed water drains through the melt water drain to a bath above the compressor (Fig. 1) and evaporates from there.

If, for some reason, the device does not defrost automatically and a layer of frost thicker than 0.5 cm forms on the rear wall, the manual defrost function can be activated: press manual defrost button, **def** will appear on display. The device compressor will stop freezing: it will run for 20 minutes. forced defrosting. After the defrost time, the device will start operating normally and the controller display will show the actual temperature inside the device.

CLEAN THE DEVICE REGULARLY.

**i** Remember that before cleaning, it is necessary to turn off with On/off button and to disconnect it from the power supply network - pull the cord plug from the electrical outlet.

- Regularly clean the hermetic gaskets of the doors. Wipe dry.
- At least once a year clean dust away from the back part of the device body and the compressor. A soft brush, electrostatic cloth or vacuum cleaner may be used for cleaning.
- REGULARLY CLEAN THE THAW WATER CHANNEL SPACE IN THE device WITH A SPECIAL CLEANER FOR THAT PURPOSE (see fig.).
- At least once a year, clean the dust accumulated on the back part of the device body, the condenser and the compressor. You can use a soft brush, electrostatic wipes or a vacuum cleaner for cleaning.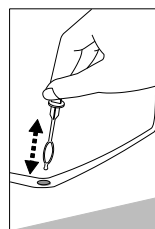

- Clean the inside, shelves and outside of the device with lukewarm water or ethyl alcohol-based cleaners or water-based household cleaners.
- Clean the glass with products intended for glass.
- DO NOT use cleaners containing abrasive particles, acid, gasoline to clean the inside and outside of the device, the door seal gasket.
- DO NOT use cloths or sponges for cleaning on rough, scrubbing surfaces.
- DO NOT use spray deicers.
- DO NOT clean with compressed air/steam.
- If the device is left empty for a long time, turn it off, defrost, clean, dry and leave the door open to prevent mold from forming inside the device.
- **Only disinfect the device with ethyl alcohol-based products.**


## HOW TO USE THE ACCESS PORT

You will need 4mm flat head screwdriver.

**ATTENTION!** Switch off the device with **On/off** button and remove the plug from the power supply socket.

1. Remove the sealing paste from the access port hole on side **A** of the compressor compartment (see Fig. 1).
2. Inside the device **B**, remove the cover from the access port hole using a screwdriver (see Fig. 2).
3. Push out the insulating material with a screwdriver (see Fig. 3).
4. Insert the sensor cable from inside the device B through the hole (see Fig. 4).
5. On the niche side A, wrap the sensor wire with insulating material and push it into the hole of the access port (see Fig. 5).
6. Additionally, seal the hole on side A with sealing paste (see Fig. 6).

After connecting the device to the mains, press and hold **On/off** button until **on** appears on screen.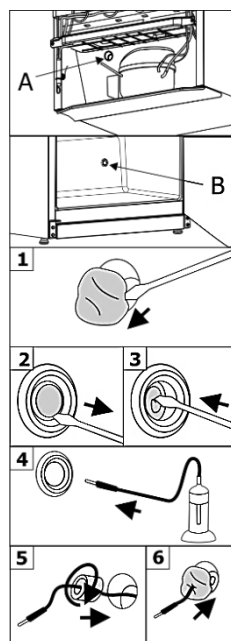

## CHANGING THE OPENING DIRECTION OF THE DOORS

We recommend changing the opening direction of the device door with an assistant. You will need: wrench 8mm and 12mm and a Phillips screwdriver.


When changing the door opening direction, it is FORBIDDEN to place the device horizontally.

**Before you start, please note the following:**


- Always ensure that the correct number of washers is used on all hinge bolts and screws.
- So that the doors can close tightly against the frame, take the utmost care that they are properly aligned during reassembly.

Do the actions below in the order they are listed:

**ATTENTION! The device is necessary to turn off with On/off button and to disconnect device from the power supply network - pull the cord plug from the electrical outlet.**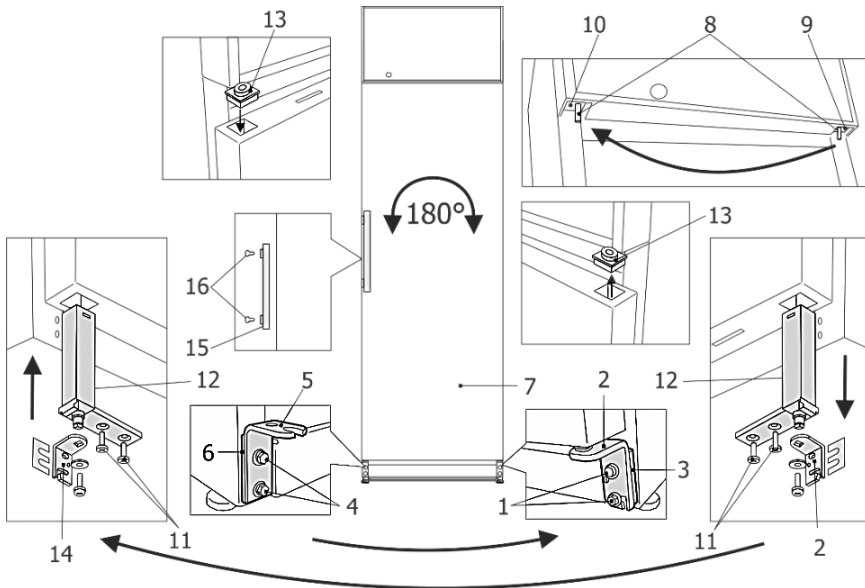

1. Turn the device at a 45° angle to the back and support it firmly.
2. Unscrew the two screws **1** and pull the door down to remove the door together with the lower bracket. and intermediate **3**.
3. Unscrew the two screws **4** and remove the bracket **5** with the spacer **6** and screw it on the opposite side of the cabinet.
4. Unscrew the axle **8** from the upper bracket **9**, (wrench 12mm) and screw into bracket **10** on the other side of the cabinet .
5. Lay the removed door on a flat surface and prepare it for left opening:
  - a) Unscrew two self-tapping screws item **11** and pull out the closing mechanism **12** from the bottom of the door.
  - b) Remove cover **13** from the top of the door and prepare a hole for inserting the door closing mechanism: remove the insulation material so that the hole for the mechanism is 100 mm deep.
  - c) Install the closing mechanism **12** and the door bushing **13** in the reverse order, changing their places.
  - d) Remove the right opening bracket **2** from the closing mechanism. and install the left opening bracket **14** in the set. Before screwing the bracket to the closing mechanism, orient the bracket so that when the door is closed there is a 15-20° opening of the door.
6. Place the door **7**, after turning it 180°, putting the pin **8** on the upper bracket **10**.

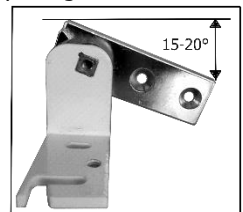7. Screw the lower bracket **14** with a set of spacers **3** on the opposite side of the cabinet.
8. Unscrew the fixing screws **16** and fix handle **15** in the holes above. Plug the loose holes with the blinds which can be found in the device packing kit bag.

If the door fits unevenly after changing the direction, you can adjust it:

- adjust the even position of the door by sliding the lower bracket left / right;
- Ensure that the door fits snugly by changing the amount of spacers **3** under the bracket **14**.

**NOTE.** Once the unit has been positioned, wait for about **30 minutes** before it is connected to the mains.

If you tilt the device at an angle of more than 45 ° when changing the direction of the door opening, allow it to stand upright for **4 hours** before connecting it to the mains.

## **OPERATION PROBLEMS AND THEIR SOLUTIONS**

- **The device is plugged in to the mains but it does not work.** Check if your house electricity supply installations are in order. Check if the plug is correctly inserted into the electricity supply socket.
- **The noise has become louder.** Check if the device is standing stable, on a level place. To make it level, regulate the front feet. Check if the device is not touching any furniture and if no part of the refrigeration system at the back part of the device body is touching the wall. Pull the device away from any furniture or walls. Check if the cause of increased noise isn't due to bottles, cans in the device that might be touching each other.
- **Water has appeared at the bottom of the compartment.** Check whether a thaw water channel space isn't blocked. Clean the thaw water channel space with a cleaner intended for that purpose.
- **When the device door is opened, the rubber sealing gasket pulls out.** The gasket is smeared with sticky foodstuffs (fat, syrup). Clean the sealing gasket and the groove for it with warm water containing soap or dishwashing detergent and wipe dry. Put the rubber sealing gasket back into the place.
- **High temperature in device, the compressor operation pauses are short.** Ascertain whether the device door/doors are shutting tightly, whether the door wasn't kept open longer than necessary when taking out or putting in foodstuffs or whether a large amount of warm food was placed in the device.
- **The top of the device has condensation on it.** The ambient relative air humidity is above 75 %. Ventilate the room where the device stands and if possible remove the cause of the humidity.
- **The temperature inside the device exceeds the set permissible limit.** Device (evaporator) does not have time to defrost automatically. Activate the manual defrost function: press the manual defrost button, the indication **DEF** will appear on the display. After the defrost time, the device will start working in normal mode and the actual temperature inside the device will be displayed on the controller display.

**Remarks on the device operating noise.** As the device operates and goes through the freezing cycle, various noises will be emitted. This is normal and is not a sign of any malfunction. As the refrigerant circulates around the refrigeration system, it causes sounds like murmuring, bubbling, or rustling. Louder sounds like popping or clicking may be heard for a short time as the refrigerator's compressor switches on.

## ENVIRONMENTAL PROTECTION INFORMATION

**Important Note:** This device is marked according to the European directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE). By ensuring this product is disposed of correctly, you will help prevent potential negative consequences for the environment and human health, which could otherwise be caused by inappropriate waste handling of this product.

The symbol on the product, or on the documents accompanying the product, indicates that this device may not be treated as household waste. Instead it shall be handed over to the applicable collection point for the recycling of electrical and electronic equipment.

For more detailed information about treatment, recovery and recycling of this product, please contact your local authority, your household waste disposal service or the retailer where you purchased the product.

Our packaging are made of environmentally friendly materials, which can be reused:

- The external packaging is made of cardboard/foil
- The FCKW free shape of foamed polystyrene (PS)
- Polyethylene (PE) foils and bags

The refrigerants and foaming agents, which are 100% free of FCKW and FKW have been used for manufacturing our product. Therefore we are helping to protect the ozone layer and prevent the increase of greenhouse effects. Also, the innovative technology and environmentally friendly insulation help in reducing energy consumption.

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AB SNAIGĖ